

Description of the Control Program CU30-0.exe


1. Connect the controller CU 30 with the piezo stages.
2. Connect the controller CU 30 with the computer via USB cable.
3. Start the program CU30-0.exe

How to start the program CU30-0.exe?

Unpack the folder "CU30-0.zip" to a folder at your computer.

Start the program "CU30-0.exe" by double-clicking.

The following screen will appear:

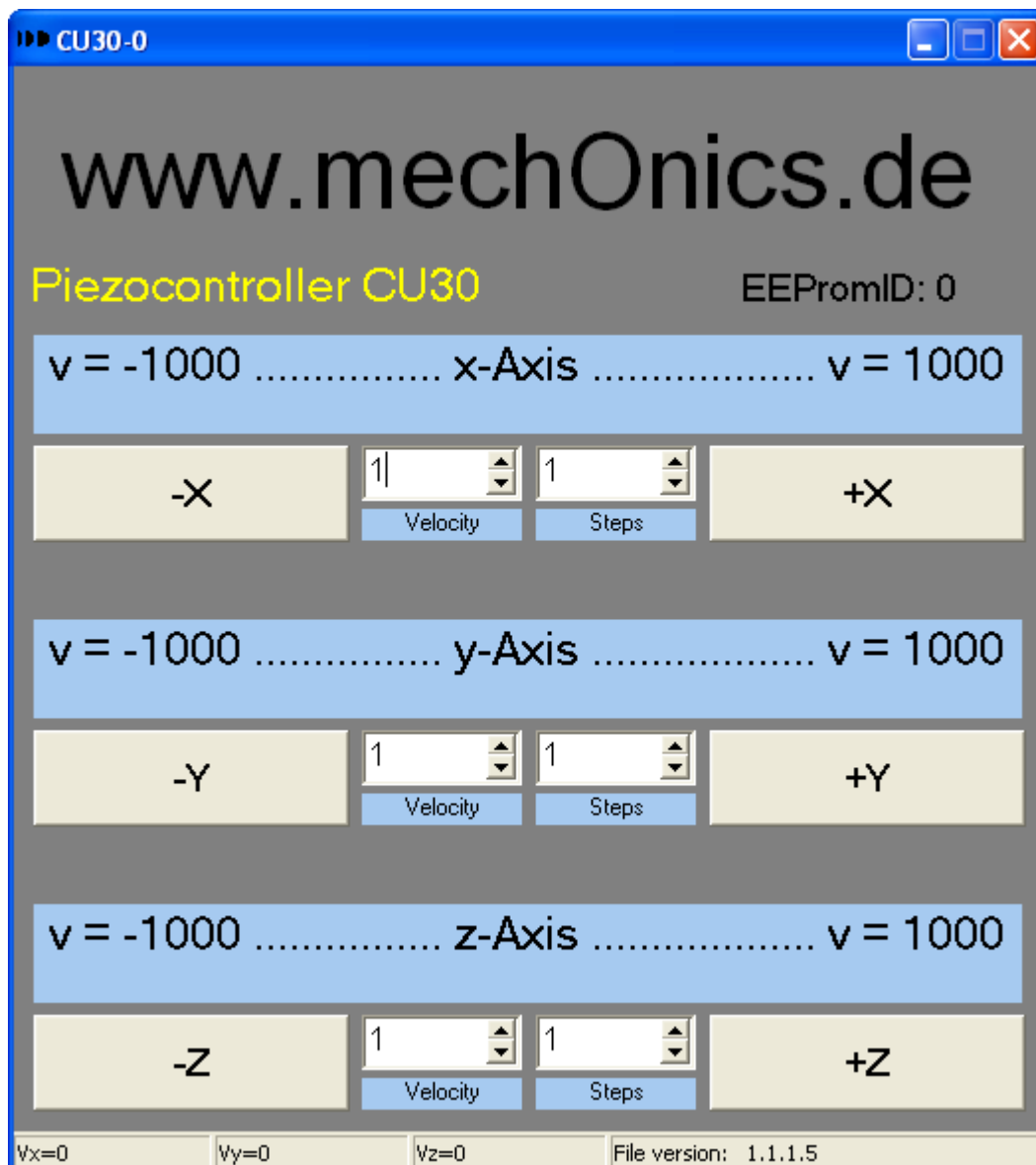

1. Swiping movement:

When you are moving with the mouse at the x, y or z - axis there will appear a double-sided arrow. By holding down the left mouse button, you can move the corresponding axis.

2. Pre-adjusted movement:

For each axis the values for velocity (1 to 1000) and steps (1 to 1,000,000) can be chosen. By clicking to the corresponding button (-X, +X, -Y, +Y, -Z or +Z) the selected number of steps will be done with the selected velocity.